

### **INSTRUCTIONS FOR SUPERFROST PLUS SLIDE USE:**

- Float your 2 to 5 micron thick tissue section on a pre-heated distilled water bath. Do not add glue or subbing solution to the bath. The Plus slide pretreatment eliminates the need for these compounds. (Note: you may try substituting tap water for distilled water. However, if you begin losing tissue sections, use only distilled water.)
- Mount your tissue sections carefully. Tissue bonding begins quickly on Superfrost Plus slides.
- Dry slides completely at room temperature by vertically draining them before heating them in an oven.