

Instructions for Superfrost Plus Slide Use

1. Float your 2 to 5 micron thick tissue section on a pre-heated distilled water bath. Do not add glue or subbing solution to the bath. The Plus slide pretreatment eliminates the need for these compounds. (Note: you may try substituting tap water for distilled water. However, if you begin losing tissue sections, use only distilled water.)
2. Mount your tissue sections carefully. Tissue bonding begins quickly on Superfrost Plus slides.
3. Dry slides completely at room temperature by vertically draining them before heating them in an oven.