

# Brain Research Laboratories

DIV. OF CAMBRIDGE INTELLIGENT SYSTEMS, INC.

P.O. BOX 88, WABAN, MASS. 02468

TEL.: (617) 965-5544

FAX: (617) 965-6220

## CHROME.-ALUM.-GELATIN COATED SLIDES.

### PREPARATION OF SLIDES;

1. Soak in 95% alcohol for at least 1/2 hr.  
Rinse twice in deionized H<sub>2</sub>O  
Rinse twice in distilled H<sub>2</sub>O
2. Dip in chrome alum solution, drain and let dry for about 1 hr.

### CHROME.-ALUM.-GELATIN SOLUTION.

2.5 gm Gelatin  
500 ml distilled H<sub>2</sub>O  
0.25gm Chrome-Alum (chromium potassium sulfate)

Use a warm water bath to help dissolve gelatin.  
Cool and add chrome-alum.  
If the solution, normally green, turns blue discard.

